QUANTOFIX® Chromat

de

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QUANTOFIX® Chromate

en

Pack content:
1 aluminum container with 100 test strips
2 test tubes
2 bottles CrO_4^{2-} -1

Measuring range:
Visually
3 – 100 mg/L CrO_4^{2-}
Reflectometrically
3 – 100 mg/L CrO_4^{2-}

Color gradation:
0 · 3 · 10 · 30 · 100 mg/L CrO_4^{2-}

Safety precautions:
Information regarding safety can be found on the box' label and in the safety data sheet. You can download the SDS from www.mn-net.com/SDS.

General indications:
Remove only as many test strips as are required. Close the container immediately after removing a strip. Do not touch the test field.

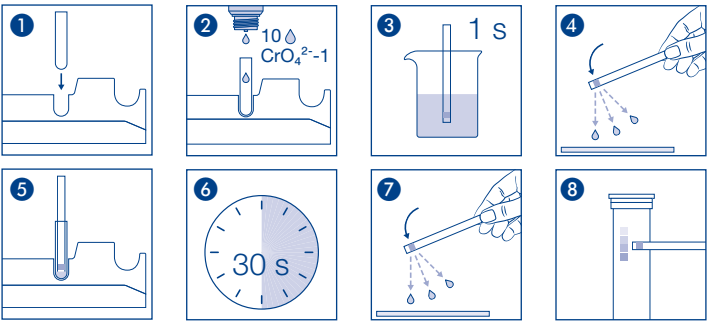
Additional indications:
If chromium is present as Cr^{3+} , it has to be converted to chromate prior to the test as follows: add to 10 mL of the test solution 1 spatula tip of potassium peroxodisulfate and 1 drop of a 0.1 mol silver nitrate solution, boil for 10 minutes and cool again to room temperature. Then, carry out the test as indicated in the instructions.
Conversion factor: $\text{Cr}^{2+} = \text{CrO}_4^{2-} \times 0.45$

- Instructions for use:**
- Place the test tube into the cavity of the thermoformed mould.
 - Add 10 drops CrO_4^{2-} -1 (nitric acid).
 - Dip the test strip into the neutral to acid test solution for 1 second.
 - Shake off excess liquid.
 - Place test strip into the filled test tube.
 - Wait 30 seconds.
 - Shake off excess liquid.
 - Compare with the color scale.
- If chromate ions are present, the test field turns violet.

Interferences:
Interferences caused by molybdates are eliminated by adding 1 spatula tip of oxalic acid to the strong acid solution. Iron(III) ions are eliminated by precipitation with 32 % sodium hydroxide solution.

Disposal:
Information regarding disposal can be found in the safety data sheet. You can download the SDS from www.mn-net.com/SDS.

Storage:
Avoid exposing the strips to sunlight and moisture. Keep container cool and dry (storage temperature 4 – 30 °C). If correctly stored, the test kit may be used until the use-by-date printed on the packaging.

MACHEREY-NAGEL GmbH & Co. KG · Valencienner Str. 11 · 52355 Düren · Germany
Tel.: +49 24 21 969-0 · info@mn-net.com · www.mn-net.com

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MACHEREY-NAGEL GmbH & Co. KG · Valencienner Str. 11 · 52355 Düren · Germania